

Detection of Human IgG Using a Luminescence-based Proximity Assay

Using Spatial Proximity Analyte Reagent Capture Luminescence (SPARCL) Assays

Paul Held Ph. D., Laboratory Manager, Applications Department, BioTek Instruments, Inc., Winooski, VT

Homogeneous proximity assays allow the quantitation of specific analytes from a complex matrix containing numerous contaminants without the need to use wash steps. Because these assays are performed in solution without binding to a solid surface, reaction kinetics are more rapid than conventional ELISA based assays. Here we describe the use the Synergy™ Neo Multi-Mode Microplate Reader in conjunction with the Spatial Proximity Analyte Reagent Capture Luminescence (SPARCL) assay technology to quantitate human IgG.

Introduction


Biomedical research requires that numerous analytes be quantitated, often in large numbers of experimental samples. Towards that end, a number of different homogeneous proximity assays have been developed to quantitate specific analytes from a complex sample matrix without the need for wash steps to remove unwanted contaminants. Eliminating the necessity to remove unbound contaminants not only saves considerable amounts of time, but also reduces the amount of instrumentation needed to automate the assays. These assays share a common feature in that signal is only generated in the presence of analyte, resulting in low background signal and excellent S/B ratios. As the name implies, this is accomplished by bringing two molecules together in close enough proximity such that an energy transfer can take place, with the analyte serving as the bridge. Typically antibodies serve as the means for analyte specificity. These features make them the preferred assay type for many screening type experiments.

A number of different homogeneous proximity assays have been developed. For example, the scintillation proximity assay (SPA) depends on the close proximity of a radio-labeled molecule to a bead containing a scintillant for signal [1]. Bioluminescence resonance energy transfer (BRET) requires that a bioluminescent moiety to be close enough to a green fluorescent protein molecule to allow the transfer of energy [2]. Time resolved fluorescence resonance energy transfer (TR-FRET) and HTRF® require that two fluorescent molecules be in close proximity, with

the donor fluor having an exceptionally long decay half-life [3-4]. Amplified Luminescent Proximity Homogeneous Assay (ALPHA) Screen assays use donor beads that generate singlet oxygen that will interact with acceptor beads located within 200 nm to generate a chemiluminescent signal [5].

There are also several nucleic acid based proximity assays. PINCER technology uses fluorescently labeled complimentary single stand nucleic acids attached to antibodies [6]. Proximity ligation assays (PLA) are based on the amplification of a DNA sequences which is dependent on the proximity of two antibodies labeled with different oligonucleotides. They form a circular DNA molecule with the help of a connector oligonucleotide, which can be detected by rolling circle amplification [7]. Proximity extension assays use target-specific antibody pairs linked to DNA strands that upon simultaneous binding to the target analyte create a real-time PCR amplicon in a proximity-dependent manner enabled by the action of a DNA polymerase [8].

Spatial Proximity Analyte Reagent Capture Luminescence (SPARCL) technology is a proximity dependent, chemiluminescent detection technique. In the assay, a chemiluminescent substrate (acridan) is brought into the proximity of an oxidative enzyme (horseradish peroxidase, HRP) through the specific antigen/antibody interaction (Figure 1). A flash of light proportional to the quantity of analyte present in the sample is generated upon addition of a trigger solution containing

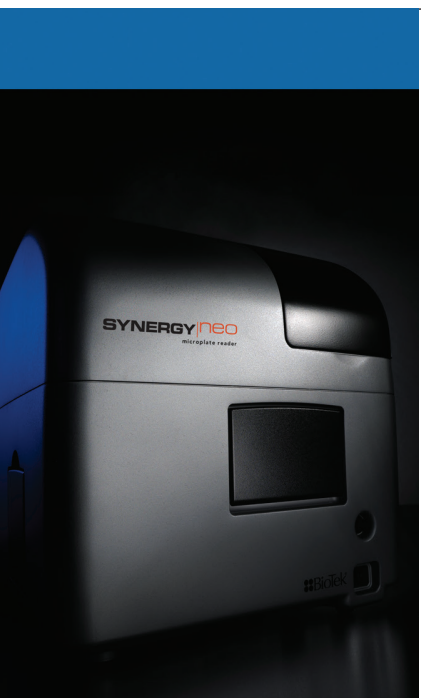

Key Words:

Luminescence

Proximity Assay

SPARCL

IgG

Homogeneous Assay

H_2O_2 and para-hydroxycinnamic acid (pHCA) enhancer. Heterocyclic compounds such as pHCA have been shown to increase light emission from HRP based luminescent reactions greater than 1000 fold [9-10]. Data from similar reactions with luminol has resulted in the proposal that enhancers serve as a conduit for the transfer of free radicals from HRP to the luminescent agent (i.e. acridan) [11]. Here we present data demonstrating SPARCL technology, a homogeneous assay based on a bridging assay format where all reagents and sample are in solution.

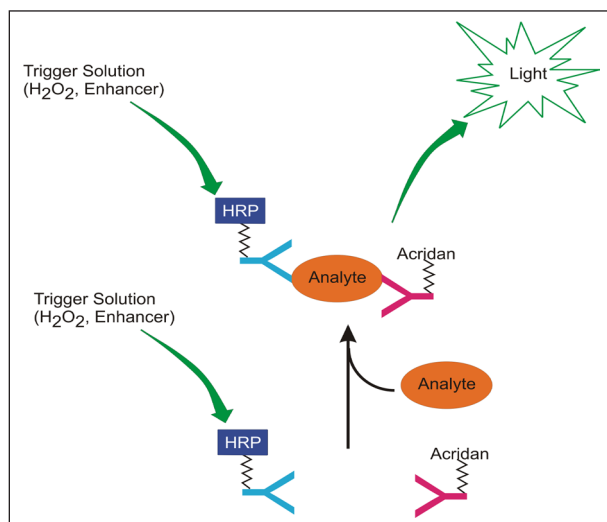


Figure 1. SPARCL reaction. The oxidative enzyme HRP is in close [proximity with the chemiluminescent substrate acridan only when both antibodies are bound to the analyte.

Horseshradish peroxidase (HRP) catalyzes the hydrolysis of hydrogen peroxide (H_2O_2) and in the process generates free radicals which oxidize substrates that are nearby. However these reactive molecules have a limited lifespan and can only diffuse short distances before they are destroyed by interacting nonspecifically with proteins or lipids. Acridan molecules in close proximity can serve as such a target for attack by the newly generated free radicals. In doing so acridans will form a dioxetone intermediate that decays to a excited acridone molecule that will emit light photons as it returns to its stable state (Figure 2). Because free radicals are highly reactive, only acridan molecules in close proximity to the source of free radicals will result in the emission of specific light photons. Only when an analyte has both a HRP conjugate and an acridan conjugate bound will the two moieties be in close enough proximity for the reaction to take place. Unbound antibodies in solution will not result in the necessary reaction proximity for a luminescent reaction to occur.

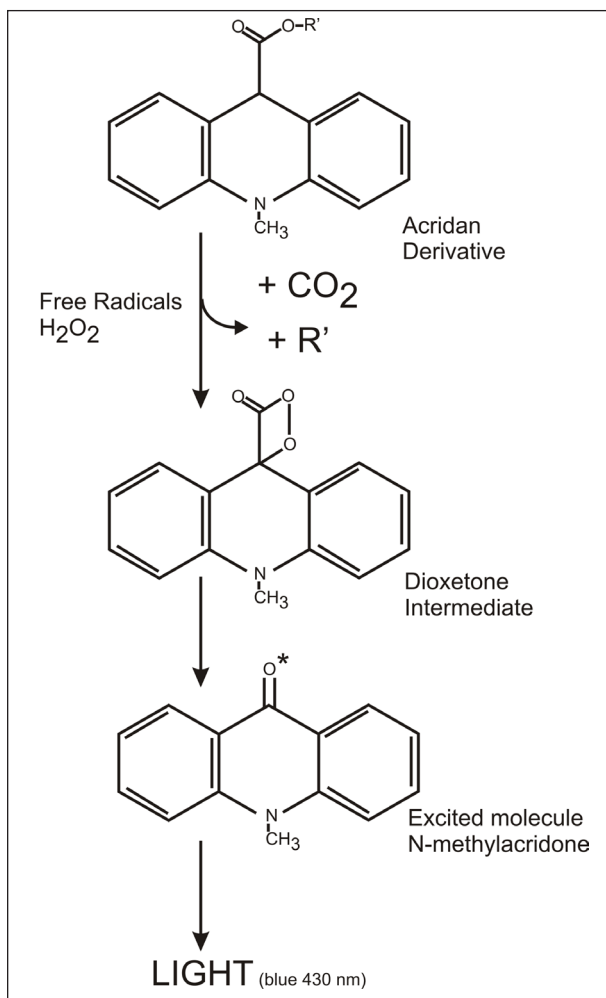


Figure 2. Acridan chemistry. An example of acridan chemistry where reaction with free radicals results in the emission of light photons.

Materials and Methods

SPARCL assay reagents including Acridan-labeled JDC-10 antibody, HRP-conjugated JDC-10 antibody, human IgG, Trigger solution, and background reducing agent were provided by Lumigen (Southfield, MI). Normal Rat serum (P/N 10710C) was obtained from Life Technologies (Carlsberg, CA) and solid white 96-well microplates (P/N 3912) and plates seals (P/N 6569) were from Corning Life Sciences (Tewksbury, MA).

SPARCL proximity assays were set up as follows. Working solutions of HRP and acridan labeled antibodies were made the day of the assay. HRP-conjugated JDC-10 antibody stock (50 $\mu\text{g}/\text{mL}$) was diluted 1:800 and acridan-labeled JDC-10 antibody stock (500 $\mu\text{g}/\text{mL}$) was diluted 1:78 with assay buffer (PBS with 0.1% BSA). Human IgG stock (10 mg/mL) was diluted to make a series of standards (0-4000 ng/mL) using normal rat serum as the diluent.

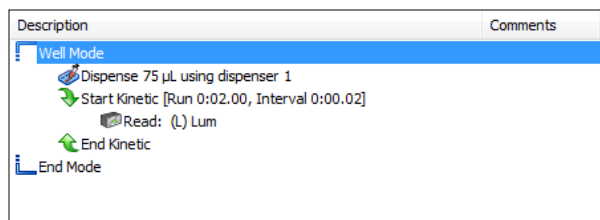


Figure 3. Dispense and Read Process. Luminescent signal is captured over 2 seconds after dispensing of trigger/enhancer reagent.

Equal amounts of HRP and acridan labeled antibody working solutions were combined and 50 μL of the mixture was aliquoted into the wells of the microplate. Samples and standards (25 μL) were then pipetted into wells of the microplate and the plate was sealed with an aluminum adhesive sealer, and incubated on a plate shaker for 30 minutes at room temperature. After incubation 4 μL of background reducing agent was added to each well and the plate was positioned on the plate reader. Luminescence was determined using a Synergy™ Neo Multi-Mode Microplate Reader (BioTek Instruments, Winooski, VT). Reactions were initiated by the injection of 75 μL of trigger reagent using the reader's reagent injectors. The reactions were measured kinetically in well mode every 0.02 seconds for a total of 2 seconds per well. Sensitivity was set to 175 and the data integral recorded (Figure 3).

Results

Kinetic luminescence data was captured using Gen5™ Data Analysis Software. When the mean signal at each time point is plotted for three different human IgG concentrations (2000, 300, 0 ng/mL) marked differences in the plots are observed. Samples containing IgG have substantial signal that diminishes over the kinetic interval, while the 0 ng/mL sample has virtually no signal (Figure 4). These data also show how rapid the luminescent reaction is.

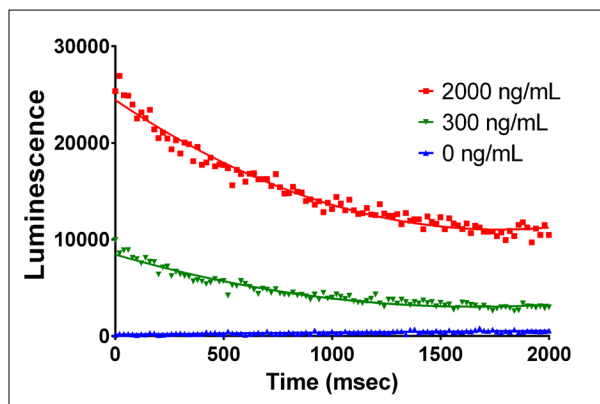


Figure 4. Kinetic luminescence of SPARCL assay for IgG at different concentrations. Kinetic data was captured every 20 msec for a period of 2 seconds. Each data point represents the mean of 10 determinations with a second polynomial regression analysis.

When the integral or area under the curve is calculated and plotted a linear relationship between analyte concentrations is observed (Figure 5). This relationship can be described using a 4-parameter logistics fit of the data, with a Z'-factor value in excess of 0.9 between the 0 and 1500 ng/mL standards.

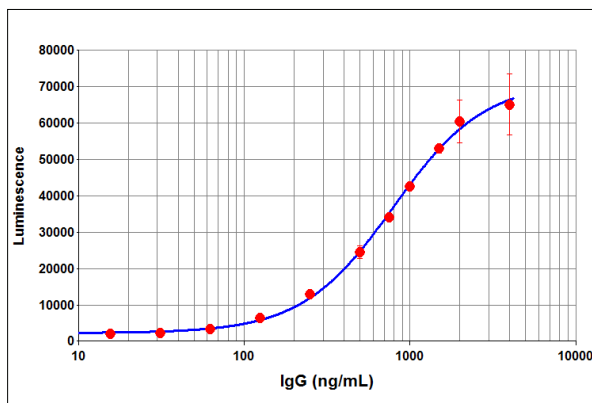


Figure 5. SPARCL IgG Titration Curve. Various concentrations of human IgG were assayed using SPARCL reagents specific to IgG. Data represents the mean and standard deviation of the integral of kinetic luminescent measurements for 4 replicates.

Discussion

These data demonstrate that the Synergy Neo is capable of making the luminescent determinations necessary for SPARCL assays. The Synergy Neo, as well as other BioTek multi-mode readers, provides a number of features that make them amenable to rapid kinetic flash-type luminescent assays. Foremost is the optional reagent dispenser module that allows for automated dispensing of two different reagents into the wells of the microplate while the plate is located in the read chamber. This feature allows for the addition of trigger reagent and enhancer and almost immediate detection of the luminescence. Gen5 Data Analysis Software (BioTek Instruments) controls reader functions, collects and stores generated data, and perform data reduction. A number of different curve fits are available to describe data or to be used as standards curve for interpolation to calculate unknown sample concentrations.

References

1. Y-W, R.T. Cummings, L. Wu, S. Zheng, P.M. Cameron, A. Woods, D.M. Zaller, A.I. Marcy, and J.D. Hermes (1999) Homogeneous Proximity Tyrosine Kinase Assays: Scintillation Proximity Assay versus Homogeneous Time-Resolved Fluorescence. *Analytical Biochemistry* **269**: 94-104.
2. Ward, W. W., Cormier, M. J., (1979). An energy transfer protein in coelenterate bioluminescence. Characterization of the Renilla green-fluorescent protein. *J. Biol. Chem.* **254**:781-788.
3. Degorce, F., A. Card, S. Soh, E. Trinquet, G. P. Knapik, and B. Xie (2009) HTRF: A Technology Tailored for Drug Discovery- A Review of Theoretical Aspects and Recent Applications, *Curr. Chem. Genomics*, **3**:22-32
4. Comley J. (2006) TR-FRET Based Assays- getting better with age *Drug Discovery World Spring* pp.22-37.
5. Eglén, R.M., T. Reisine, P. Roby, N. Rouleau, C. Illy, R. Bossé, and M. Bielefeld (2008) The Use of AlphaScreen Technology in HTS: Current Status, *Curr. Chem. Genomics*, **1**:2-10.
6. Heyduk, E. B. Dummit, Y-W. Chang, and T. Heyduk (2008) Molecular Pincers-new Antibody-based Homogeneous Protein Sensors, *Anal. Chem.* **80**:5152-5159.
7. Fredriksson S, M. Gullberg, J. Jarvius, C. Olsson, K. Pietras, S.M. Gustafsdottir, A. Ostman, and U. Landegren (2002) Protein detection using proximity-dependent DNA ligation assays. *Nat. Biotechnol.* **20**:473-477.
8. Lundberg, M., A. Eriksson, B. Tran, E. Assarsson and S. Fredriksson (2011) Homogeneous antibody-based proximity extension assays provide sensitive and specific detection of low-abundant proteins in human blood. *Nucleic Acids Research*, p1-8. doi:10.1093/nar/gkr424
9. García-Sánchez, F. A. Navas Díaz, and J.A. González García (1997) Relation between the structure of some heterocyclic derivatives and other compounds, and their effects as enhancers or inhibitors of the luminal-H₂O₂-horseradich peroxidase chemiluminescence, *J. of Photochemistry and Photobiology A: Chemistry*, **105**:11-14.
10. Thorpe, G.H.G., L.J. Kricks, S.B. Mosely, T.P. Whitehead (1985) Phenols as Enhancers of the Chemiluminescent Horseradish Peroxidase-Luminol-Hydrogen Peroxide Reaction: Application in Luminescence-Monitored Enzyme Immunoassays, *Clinical Chemistry*, **31/8**:1335-1341.
11. Garcia Sánchez, F., A. Díaz Diaz, and J.A. González García (1995) P-phenol Derivatives as Enhancers of the Chemiluminescent Luminol-Horseradish Peroxidase-H₂O₂ Reaction: Substituent Effects, *J. Luminescence*, **65**:33-39.